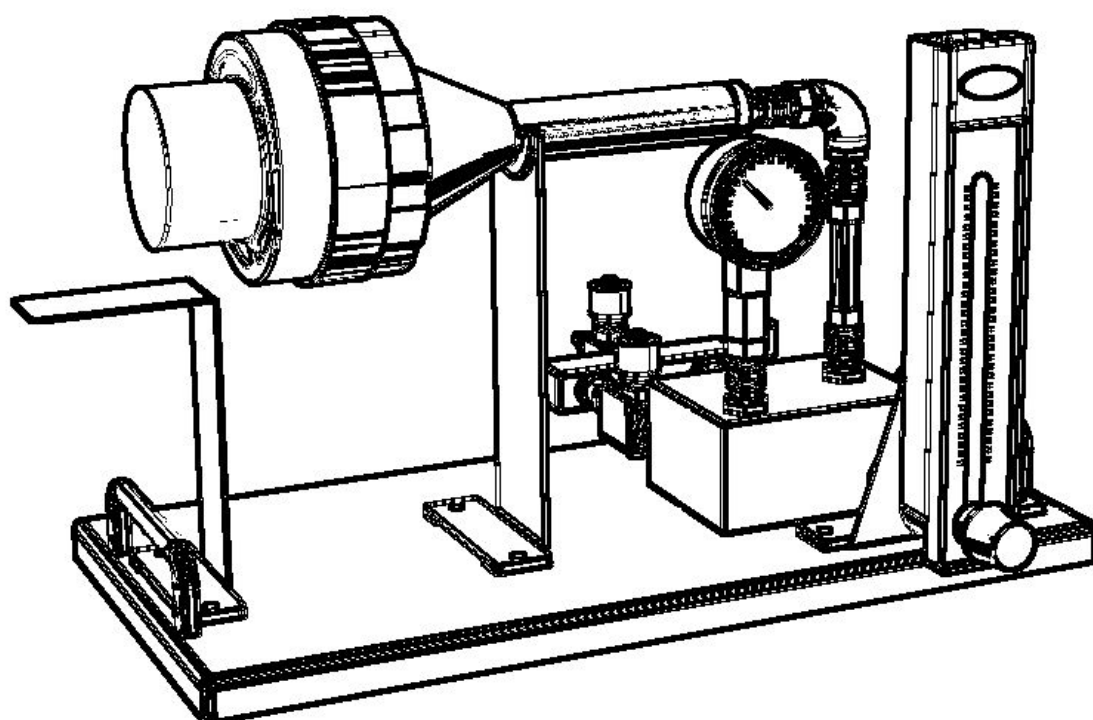
VWR® Pinocchio Super

INSTRUCTION MANUAL

European catalogue numbers:

710-0950	Pinocchio Super Petri
710-0949	Pinocchio Super Contact
710-0955	Pinocchio Super Petri Validated
710-0951	Pinocchio Super Contact Validated

Version: 2.2
Issued: July 2016

Legal address of manufacturer

Europe

VWR International S.r.l.

Via San Giusto, 85

20153 Milano (MI)

Italy

 + 39 023320311

<http://it.vwr.com>

Country of origin: Italy

Table of contents

Warning	4
Safety information	4
Installation	5
Intended use	7
Specifications	7
Instructions for	8
Maintenance	10
Technical service	11
Warranty	11
Equipment disposal	12

Warning

PLEASE READ THE INFORMATION CONTAINED IN THIS MANUAL BEFORE USING THE UNIT.
THE MANUFACTURER DOES NOT ACKNOWLEDGE ANY RESPONSIBILITY FOR AN IMPROPER USE
OF THIS EQUIPMENT AND NOT FOLLOWING THESE DIRECTIONS FOR USE.

Using this manual will help you use this instrument efficiently. The manual should be kept in a convenient place near the instrument for ease of use.

Safety information

Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

- Use the device just for the indicated purposes
- The device must be correctly installed according to manual instructions, before starting any operation
- Use original spare parts and accessories for any replacement
- If this device is used with extremely high pressurised gas you must connect a reducer to the inlet in order to have a maximum of 6 bar to the inlet

Avoiding damage to the instrument

Be aware that pressure applied to the inlet is lower than 6 bar.

Avoiding damage to operator

Do not use this device with gas with an oxygen concentration more than 21%.

Do not use this device with gas harmful for humans or the environment.

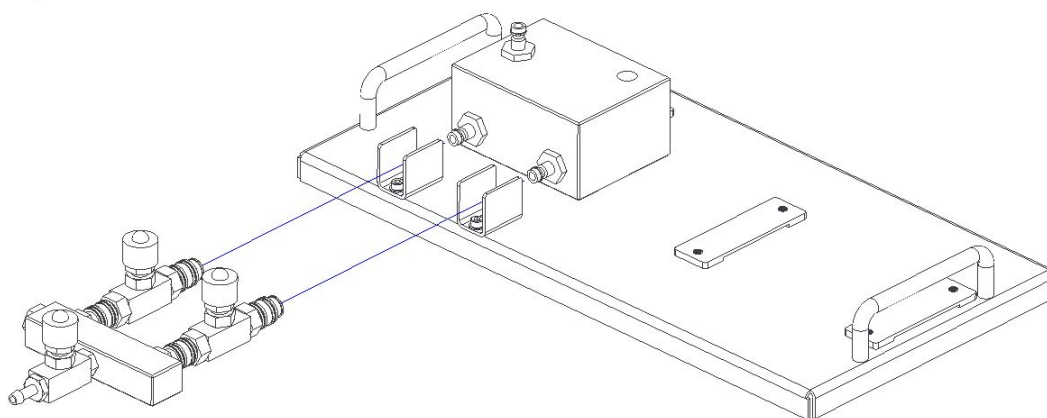
Installation

The compactness of “Pinocchio Super II” sampler enables the device to be located where space is limited. The simple construction coupled with simple operation enable accurate reports on air–gas microbiology quality. The unit can be completely dismantled for easy cleaning and autoclave sterilisation. The pressure gauge and the airflow meter are not autoclavable. Follow the schematic drawing for assembly. All connections are “quick disconnect”.


Connect the unit to the desired point of sampling using suitable sterile tubing. The inlet gas valve should be in a closed position.

ASSEMBLING


- 1) Connect the inlet splitter to the basement according to the following picture.

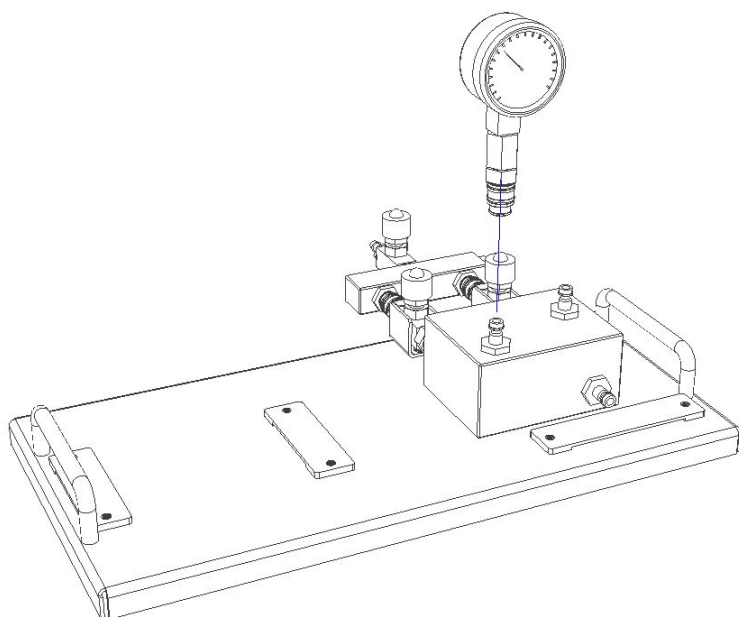
- 2) Connect the pressure gauge to the base according to the following picture.

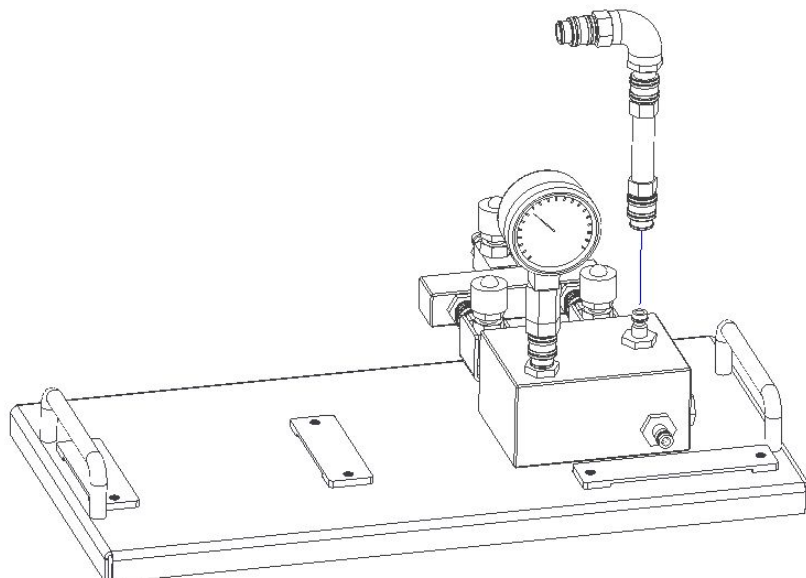
3) Connect the elbow arm to the base according to the following picture.

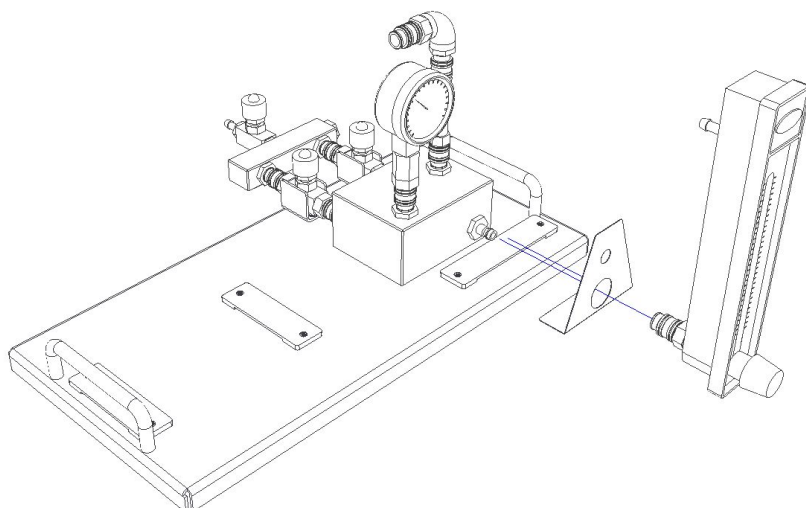4) Connect the flow meter and its holder to the base according to the following picture.

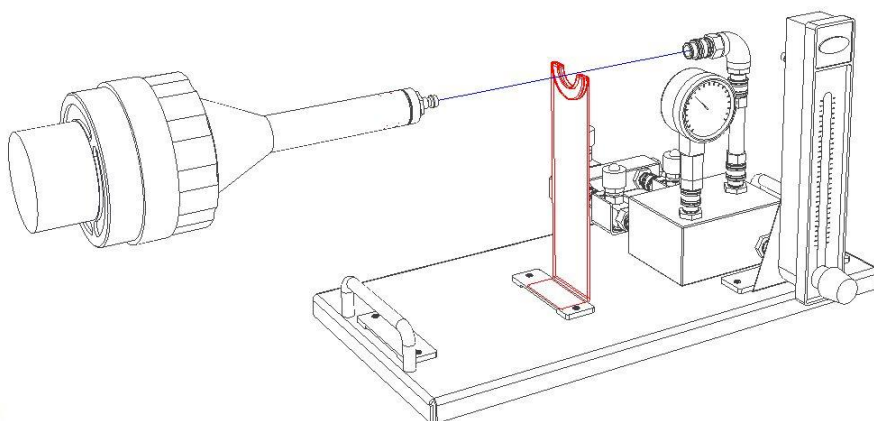5) Connect the sampling arm and its holder to the base according to the following picture.

Intended use

The “Pinocchio Super II” microbiological gas sampler is a portable instrument for determining the level of existent microbiological contaminants in the compressed gas of aseptic manufacturing operations.

For laboratory use only. Not intended for any animal or human therapeutic or diagnostic use.

Specifications

HxWxD	245x400x200 mm
Weight	5 kg
Construction	Stainless steel / anodized aluminium
Airflow meter	50/400 litres
Operating temperature	10 40 °C
Electrical power supply	N/A
Compressed air	Max. inlet pressure 6 bar

Overview

The Pinocchio Super II microbiological gas sampler is a portable instrument for determining the level of existent microbiological contaminants in the compressed gas of aseptic manufacturing operations.

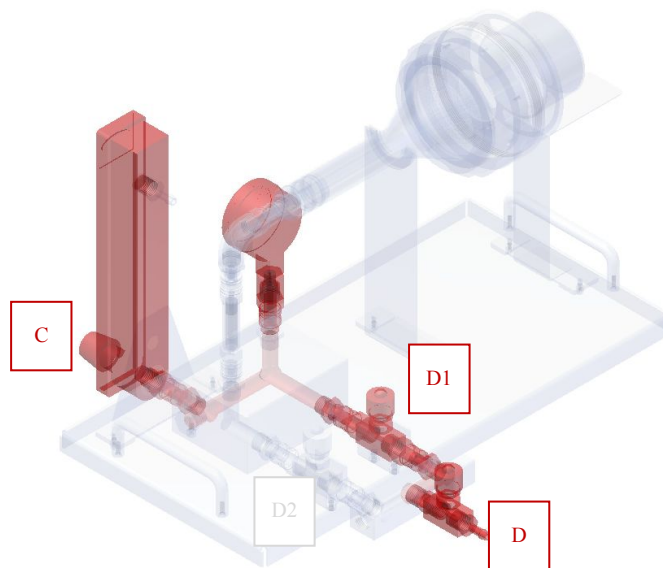
An official validation of the Pinocchio Super II air sampler is required as part of an inspection. The flow meter should be calibrated by an official agency at regular intervals (usually every year). Official validation of the flow meter can also be performed by VWR INTERNATIONAL customer support service.

Instructions for use

PREPARATION


- 1) Connect the unit to the desired point of sampling using suitable sterile tubing. The inlet gas valve should be in a closed position.
- 2) Insert an identified contact plate or Petri dish with sterile medium in the Pinocchio funnel housing and screw it to the conical head. Do not connect to the Pinocchio Super II.

AIRFLOW ADJUSTMENT

- 3) Fully open the flow meter knob (pos. "C" Figure 1).
- 4) Fully close the D and D2 valves.
- 5) Fully open the D1 valve.
- 6) Open the line of compressed air or gas.
- 7) Start opening the D valve slightly, ensuring that the manometer is not indicating a pressure greater than 6 bar.
- 8) Open the D valve until you see the desired airflow on the flow meter, the suggest value is 100 l/min (210 SCFH). The purpose of this operation is to obtain a known volume of air, without contaminating the sampling circuit.

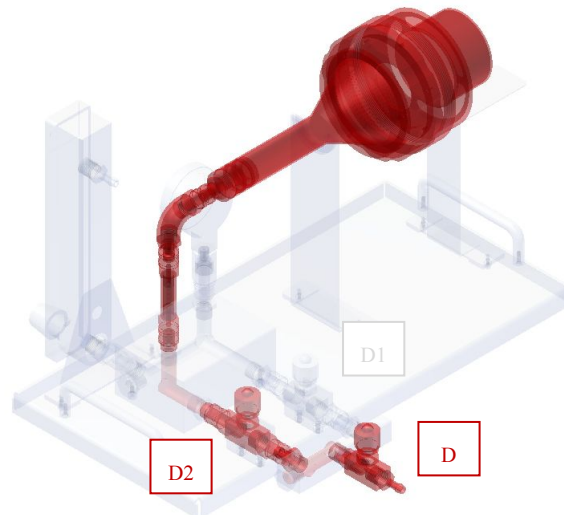COMPRESSED GAS SAMPLING

- 9) Connect the Pinocchio funnel to the Pinocchio Super II.
- 10) Fully close the (D1) valve.
- 11) Fully open the (D2) valve and start the count down.
- 12) Close the inlet gas valve after 10 minutes (or when 1000 litres of air is collected).
- 13) Unscrew the head and transfer the contact plate (RODAC) or Petri dish to a laboratory for incubation.

The Pinocchio Super II is now ready for a new test.

Count the Colony Forming Units in the plate at the end of incubation time and report the results as CFU/1000 litres of gas/compressed air.

SAMPLING DATA RECORDING

If the sampling data needs to be recorded, replace the plate housing with a SAS Super air sampler and start from point 4 to 6. The time will be regulated by the sampler timer.

CALIBRATION

The airflow meter can be officially calibrated each year.

CLEANING & STERILISATION

The unit can be dismantled for easy cleaning and autoclave sterilisation.

IMPORTANT NOTE!

AIRFLOW METER AND PRESSURE GAUGE ARE NOT AUTOCLAVABLE.

CONVERSION TABLE “SCFH” TO LITRES PER MINUTE

The meters airflow readings should be translated into litres per minute. Follow the reported table.

SCFH	l/min
150	70,8
160	75,5
170	80,2
180	85,0
190	89,7
200	94,4
210	99,1
220	103,8
230	108,5
240	113,3
250	118,0
260	122,7

Maintenance

Daily cleaning of the instrument will prevent most future faults. It is possible to clean the instrument with a mild detergent and a soft wet cloth.

WARNING

Do not use any cleaning product containing halogens or other aggressive substance to clean the instrument. Usage of these products can cause rust.

Technical service

Web resources

Visit the VWR website at vwr.com for:

- Complete technical service contact information
- Access to the VWR online catalogue and information about accessories and related products
- Additional product information and special offers

Contact us For information or technical assistance contact your local VWR representative or visit vwr.com.

Warranty

VWR International warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorisations necessary to run or use the product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorisation, unless any refusal is due to a defect of the product.

Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle end by handing it over to an authorised facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

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Obrigado

Local VWR offices in Europe and Asia Pacific

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India

VWR Lab Products Private Limited
135/12, Brigade Towers, 2nd Floor
Front wing, Brigade Road,
Bengaluru, India – 560 025
Tel: +91-2522-647911/922 (Mumbai)
Tel: +91-80-41117125/26 (Bangalore)
Fax +91-80-41117120
E-mail: vwr_india@vwr.com

Ireland / Northern Ireland

VWR International Ltd / VWR International
(Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: 01 88 22 222
Fax: 01 88 22 333
E-mail sales@ie.vwr.com

Italy

VWR International S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: 02-3320311/02-487791
Fax: 02-332031307/02-40090010
E-mail: info@it.vwr.com
info@it.vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: 020 4808 400
Fax: 020 4808 480
E-mail: info@nl.vwr.com

New Zealand

VWR International LP
241 Bush Road
Albany
0632 New Zealand

Norway

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: 02290
Fax: 815 00 940
E-mail: info@no.vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: 058 32 38 210
Fax. 058 32 38 205
E-mail: labart@pl.vwr.com

Portugal

VWR International - Material de
Laboratório, Lda
Edifício Neopark
Av. Tomás Ribeiro, 43- 3 D
2790-221 Carnaxide
Tel.: 21 3600 770
Fax: 21 3600 798/9
E-mail: info@pt.vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel: +65 6505 0760
Fax: +65 6264 3780
E-mail: sales@sg.vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: 902 222 897
Fax: 902 430 657
E-mail: info@es.vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: 08 621 34 00
Fax: 08 621 34 66
E-mail: info@se.vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: 044 745 13 13
Fax: 044 745 13 10
E-mail: info@ch.vwr.com

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard
Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: 0800 22 33 44
Fax: 01455 55 85 86
E-mail: uksales@uk.vwr.com