



## Circulating Baths with MX Temperature Controller

### INSTRUCTION MANUAL



#### North American Catalog Number

|   | <u>120 V</u> | <u>240 V</u> |
|---|--------------|--------------|
| MX07R-20 Refrigerating / Heating Bath   | 89171-252    | 89171-254    |
| MX7LR-20 Refrigerating / Heating Bath   | 89171-244    | 89171-246    |
| MX06S135 Stainless Steel Open Tank Bath | 89171-196    | 89171-198    |
| MX10S135 Stainless Steel Open Tank Bath | 89171-200    | 89171-202    |
| MX20S135 Stainless Steel Open Tank Bath | 89171-204    | 89171-206    |
| MX11P100 Polycarbonate Open Tank Bath   | 89171-176    | 89171-178    |
| MX17P100 Polycarbonate Open Tank Bath   | 89171-180    | 89171-182    |
| MX28P100 Polycarbonate Open Tank Bath   | 89171-184    | 89171-186    |
| MX29VB3S Viscosity Bath                 | 89171-192    | 89171-194    |
| MX29VB5R Viscosity Bath                 | 89171-188    | 89171-190    |
| MX28C135 Coliform Bath                  | 89171-208    | 89171-210    |

#### European Catalog Number

|   | <u>240 V</u> |
|---|--------------|
| MX07R-20 Refrigerating / Heating Bath   | 462-0225     |
| MX7LR-20 Refrigerating / Heating Bath   | 462-0223     |
| MX06S135 Stainless Steel Open Tank Bath | 462-0211     |
| MX10S135 Stainless Steel Open Tank Bath | 462-0212     |
| MX20S135 Stainless Steel Open Tank Bath | 462-0213     |
| MX11P100 Polycarbonate Open Tank Bath   | 462-0206     |
| MX17P100 Polycarbonate Open Tank Bath   | 462-0207     |
| MX28P100 Polycarbonate Open Tank Bath   | 462-0208     |
| MX29VB3S Viscosity Bath                 | 462-0210     |
| MX29VB5R Viscosity Bath                 | 462-0209     |
| MX28C135 Coliform Bath                  | 462-0214     |

Version: 1  
Issued: 17 May 2011



## Legal Address of Manufacturer

### **United States**

VWR International, LLC  
Radnor Corporate Center, Building One, Suite 200  
P.O. Box 6660, 100 Matsonford Road  
Radnor, PA 19087 U.S.A.  
800-932-5000  
<http://www.vwr.com>

### **Europe**

VWR International bvba  
Researchpark Haasrode 2020  
Geldenaaksebaan 464  
B-3001 Leuven  
+ 32 16 385011  
<http://be.vwr.com>

## Country of Origin

United States of America

# Table of Contents

|  |           |
|--|-----------|
| <b>Introduction .....</b>  | <b>2</b>  |
| VWR Circulating Baths with the MX Temperature Controller .....                   | 2         |
| General Safety Information .....   | 3         |
| Safety Recommendations .....   | 4         |
| Regulatory Compliance and Testing .....  | 5         |
| Unpacking Your Circulator .....  | 5         |
| Package Contents .....   | 6         |
| Controls and Components .....  | 7         |
| <b>Quick-Start .....</b>   | <b>10</b> |
| <b>Installation and Startup .....</b>  | <b>12</b> |
| General Site Requirements .....  | 12        |
| Adding Liquid to the Bath Reservoir .....  | 12        |
| External Closed Loop Circulation .....   | 13        |
| Refrigeration Control Connections (Refrigerating/Heating Circulators only) ..... | 14        |
| Electrical Power .....   | 15        |
| Controller Setup .....   | 16        |
| <b>Normal Operation .....</b>  | <b>17</b> |
| Keys and Controls .....  | 17        |
| Turning Your Circulator ON .....   | 17        |
| Main Operational Display .....   | 17        |
| Set-Up Sub-Menus .....   | 18        |
| Adjusting the Temperature Set Point .....  | 18        |
| Selecting the Temperature Unit .....   | 19        |
| Setting the Offset Calibration .....   | 20        |
| Setting the Low Limit Temperature .....  | 21        |
| Setting the High Limit Temperature .....   | 21        |
| Adjusting the Auto Cool Temperature .....  | 22        |
| Resetting the Factory Default Values .....                                       | 23        |
| Loss of Power Restart .....  | 23        |
| Tap Water Cooling .....  | 24        |
| Reservoir Cover Storage .....  | 24        |
| <b>Display Messages and Alarms .....</b>   | <b>25</b> |
| <b>Routine Maintenance and Troubleshooting .....</b>                             | <b>26</b> |
| Maintaining Clear Bath Water .....   | 26        |
| Draining the Bath Reservoir .....  | 27        |
| Checking the Over-Temperature Safety System .....                                | 28        |
| Cleaning Your Circulator .....   | 28        |
| Temperature Controller .....   | 28        |
| Bath Reservoir .....   | 28        |
| Condenser, Air Vents, and Reusable Filter (Refrigerated Models only) .....       | 29        |
| Temperature Controller Removal and Re-Installation .....                         | 30        |
| Troubleshooting Chart .....  | 32        |
| <b>Technical Information .....</b>   | <b>34</b> |
| Performance Specifications .....   | 34        |
| Reservoir Fluids .....   | 35        |
| Application Notes .....  | 36        |
| <b>Equipment Disposal (WEEE Directive) .....</b>                                 | <b>37</b> |
| <b>Technical Service .....</b>   | <b>37</b> |
| <b>Warranty .....</b>  | <b>37</b> |
| <b>Local VWR Offices in Europe and Asia Pacific .....</b>                        | <b>38</b> |

## Introduction

Thank you for choosing a VWR Circulating Bath with MX Temperature Controller. It is intended for the precise temperature control of suitable non-flammable liquids (per DIN 12876-1) in a reservoir.



**WARNING:** VWR Circulating Baths are not intended for directly controlling the temperature of foods, pharmaceuticals, medicines, or other objects which may be ingested by or injected in humans or animals. Any such objects must be isolated from contact with the bath fluid and bath surfaces.

It will take you very little time to get your new Circulating Bath installed and running. This Instruction Manual is designed to guide you quickly through the process. We recommend that you read it thoroughly before you begin.









### VWR Circulating Baths with the MX Temperature Controller

| Model Type                              | Reservoir Capacity | Temperature Range                   |                                      |
|---|--------------------|-------------------------------------|--------------------------------------|
|   |                    | °C                                  | °F                                   |
| MX07R-20 Refrigerating / Heating Bath   | 7 liters           | -20° to 135°C                       | -4° to 275°F                         |
| MX7LR-20 Refrigerating / Heating Bath   | 7 liters           | -20° to 135°C                       | -4° to 275°F                         |
| MX06S135 Stainless Steel Open Tank Bath | 6 liters           | Ambient +10° to 135°C               | Ambient +20° to 275°F                |
| MX10S135 Stainless Steel Open Tank Bath | 10 liters          | Ambient +10° to 135°C               | Ambient +20° to 275°F                |
| MX20S135 Stainless Steel Open Tank Bath | 20 liters          | Ambient +10° to 135°C               | Ambient +20° to 275°F                |
| MX11P100 Polycarbonate Open Tank Bath   | 11 liters          | Ambient +10° to 85°C <sup>(1)</sup> | Ambient +20° to 185°F <sup>(1)</sup> |
| MX17P100 Polycarbonate Open Tank Bath   | 17 liters          | Ambient +10° to 85°C <sup>(1)</sup> | Ambient +20° to 185°F <sup>(1)</sup> |
| MX28P100 Polycarbonate Open Tank Bath   | 28 liters          | Ambient +10° to 85°C <sup>(1)</sup> | Ambient +20° to 185°F <sup>(1)</sup> |
| MX29VB3S Polycarbonate Viscosity Bath   | 27 liters          | Ambient +10° to 85°C <sup>(1)</sup> | Ambient +20° to 185°F <sup>(1)</sup> |
| MX29VB5R Polycarbonate Viscosity Bath   | 17 liters          | Ambient +10° to 85°C <sup>(1)</sup> | Ambient +20° to 185°F <sup>(1)</sup> |
| MX28C135 Coliform Bath                  | 28 liters          | Ambient +10° to 135°C               | Ambient +20° to 275°F                |

1. Maximum operating temperature for polycarbonate tank; MX Controller is capable of higher temperatures.

## General Safety Information






When installed, operated, and maintained according to the directions in this manual and common safety procedures, your Circulating Bath should provide safe and reliable temperature control. Please ensure that all individuals involved in the installation, operation, or maintenance of this Circulating Bath read this manual thoroughly prior to working with the unit.

|   |  |
|---|--|
|    | This symbol alerts you to a wide range of potential dangers.   |
|    | This symbol advises danger from electricity or electric shock.   |
|    | This symbol indicates that a hot surface may be present.   |
|    | This symbol marks information that is particularly important.  |
|   | This symbol indicates alternating current.   |
|  | These symbols on the Power Switch / Circuit Breaker indicate that they place the main power supply ON / OFF.                                   |
|  | This symbol on the Power Key indicates that it places the unit in a standby mode. It DOES NOT fully disconnect the unit from the power supply. |
|  | This symbol indicates a protective conductor terminal.   |

**Read all instructions pertaining to safety, set-up, operation, and maintenance.  
Proper operation is the user's responsibility.**

## Safety Recommendations

To prevent injury to personnel and/or damage to property, always follow your workplace's safety procedures when operating this equipment. You should also comply with the following safety recommendations:

|   |   |
|---|---|
|    | <p><b>WARNING:</b></p> <ul style="list-style-type: none"><li>• This Circulating Bath is suitable only for use with Class I non-flammable fluids (per DIN 12876-1).</li><li>• Be aware of the chemical hazards that may be associated with the bath fluid used. Observe all safety warnings for the fluids used as well as those contained in the material safety data sheet.</li><li>• Use only recommended bath fluids.</li><li>• Use only non-acid bath fluids.</li></ul>   |
|    | <p><b>WARNING:</b></p> <ul style="list-style-type: none"><li>• Always connect the power cord on this Circulator to a grounded (3-prong) power outlet. Make certain that the outlet is the same voltage and frequency as your unit.</li><li>• Never operate the Circulator with a damaged power cord.</li><li>• Always turn the Circulator OFF and disconnect mains power before performing any maintenance or service.</li></ul>  |
|   | <p><b>WARNING:</b></p> <ul style="list-style-type: none"><li>• Never operate the Circulator without bath fluid in the reservoir. Periodically check the reservoir to ensure that the liquid depth is within acceptable levels. Always refill the reservoir using the same bath fluid that is already in the reservoir. Bath oil must not contain any water contaminants and should be preheated to the actual bath temperature before adding as there is an explosion hazard at high temperatures.</li><li>• Always drain all fluid from the reservoir before moving or lifting your Circulator. Be sure to follow your organization's procedures and practices regarding the safe lifting and relocation of heavy objects.</li></ul> |
|  | <p><b>WARNING:</b></p> <ul style="list-style-type: none"><li>• Always allow the bath fluid to cool to ambient temperature before draining.</li><li>• The reservoir cover, top deck, and/or external pump connections may become hot with continuous use. Exercise caution when touching these parts.</li><li>• Always keep within the 85°C maximum operating temperature limit if using a polycarbonate open bath tank.</li></ul>   |
|  | <p><b>WARNING:</b> It is the user's responsibility to properly decontaminate the unit in the event hazardous materials are spilled on exterior or interior surfaces. Consult manufacturer if there is any doubt regarding the compatibility of decontamination or cleaning agents.</p>  |

## Regulatory Compliance and Testing

This equipment is compliant with the European Directive 2002/95/EC and its latest amendments on Restrictions on Hazardous Substances (RoHS) and below the given limits of hazardous substances.

### ETL Intertek (60 Hz units)

UL 61010-1 / CSA C22.2 No. 61010-1 — Safety Requirements for Measurement, Control, and Laboratory Use; Part 1: General Requirements

UL 61010A-2-010 / CSA C22.2 No. 61010-2-010:04 — Safety Requirements for Measurement, Control, and Laboratory Use; Part 2: Particular Requirements for Laboratory Equipment for the Heating of Materials

UL 508C — Standards for Safety for Power Conversion Equipment

UL 873 — Standard for Safety for Temperature-Indicating and –Regulating Equipment

UL 427 — Standard for Safety for Refrigerating Units

CSA C22.2 No. 120-M91 — Standard for Safety for Refrigeration Equipment

### CE (50 Hz units)

EC Low Voltage Directive 2006/95/EC

EC Electromagnetic Compatibility Directive 2004/108/EC

IEC 61010-1-2001

IEC 61010-2-2001

IEC 61326:2005 / EN 61326 : 2006

## Unpacking Your Circulator

Your Circulator is packed in a special carton or cartons. You should keep the packaging, along with all packing materials, until the unit has been installed and you are certain it is working properly.



**CAUTION:** Remove any loose packing material that may have fallen into the heater/pump housing during shipping. Before powering up, check that nothing remains around the heater or Circulator pump.

We recommend that you begin using your Circulator immediately to confirm proper operation, since beyond one week you may be eligible for warranty repair only (rather than replacement). You'll find complete warranty information in the back of this manual.

In the unlikely event that the unit was damaged or does not operate properly, immediately contact the transportation company, file a damage claim, and contact the company where your Circulator was purchased.

## Package Contents

The items included with your Circulator will vary depending on which model Circulating Bath you purchased.

### Models with North American power cord:

|   | Refrigerating / Heating Circulators | Stainless Steel Open Tank Circulators | Polycarbonate Open Tank Circulators | Viscosity Baths  | Coliform Bath  |
|---|-------------------------------------|---------------------------------------|-------------------------------------|--|----------------|
| Resource Disk (with Instruction Manual) | 1                                   | 1                                     | 1                                   | 1  | 1              |
| Reservoir Lid / Bath Cover              | 1                                   | 1 <sup>1</sup>                        | 1 <sup>1</sup>                      | 5<br>(models with round opening)<br>3<br>(models with square openings) | 1 <sup>1</sup> |
| 3-ft / 0.91 m IEC to IEC Power Cord     | 1                                   | N/A                                   | N/A                                 | N/A  | N/A            |
| 6-ft / 1.82 m IEC to Mains Power Cord   | 1                                   | 1                                     | 1                                   | 1  | 1              |
| Refrigeration Control Cable             | 1                                   | N/A                                   | N/A                                 | N/A  | N/A            |
| Fitting kit for external applications   | 1                                   | 1                                     | 1                                   | N/A  | N/A            |
| Cooling coil                            | N/A                                 |                                       |                                     | 1  |                |
| Qualification Test Results              | 1                                   | 1                                     | 1                                   | 1  | 1              |
| Quick-Start Guide                       | 1                                   | 1                                     | 1                                   | 1  | 1              |

### Models with European power cords:


|  | Refrigerating / Heating Circulators | Stainless Steel Open Tank Circulators | Polycarbonate Open Tank Circulators | Viscosity Baths  | Coliform Bath  |
|--|-------------------------------------|---------------------------------------|-------------------------------------|--|----------------|
| Resource Disk (with Instruction Manual)            | 1                                   | 1                                     | 1                                   | 1  | 1              |
| Reservoir Lid / Bath Cover                         | 1<br>(2 on 45L model)               | 1 <sup>1</sup>                        | 1 <sup>1</sup>                      | 5<br>(models with round opening)<br>3<br>(models with square openings) | 1 <sup>1</sup> |
| 3-ft / 0.91 m IEC to IEC Power Cord                | 1                                   | N/A                                   | N/A                                 | N/A  | N/A            |
| 6-ft / 1.82 m IEC to Mains Power Cord              | 1                                   | 1                                     | 1                                   | 1  | 1              |
| 6-ft / 1.82 m IEC to Mains Power Cord – EU plug    | 1                                   | 1                                     | 1                                   | 1  | 1              |
| 6-ft / 1.82 m IEC to Mains Power Cord – UK plug    | 1                                   | 1                                     | 1                                   | 1  | 1              |
| 6-ft / 1.82 m IEC to Mains Power Cord – Swiss plug | 1                                   | 1                                     | 1                                   | 1  | 1              |
| Refrigeration Control Cable                        | 1                                   | N/A                                   | N/A                                 | N/A  | N/A            |
| Fitting kit for external applications              | 1                                   | 1                                     | 1                                   | N/A  | N/A            |
| Qualification Test Results                         | 1                                   | 1                                     | 1                                   | 1  | 1              |
| Quick-Start Guide                                  | 1                                   | 1                                     | 1                                   | 1  | 1              |

1. 10 and 20 liter Stainless Steel Open Tank Baths. 17 and 28 liter Polycarbonate Open Tank Baths, and the Coliform Bath also have a lidded opening that accommodates an optional cooling coil.



## Controls and Components

### MX Controller

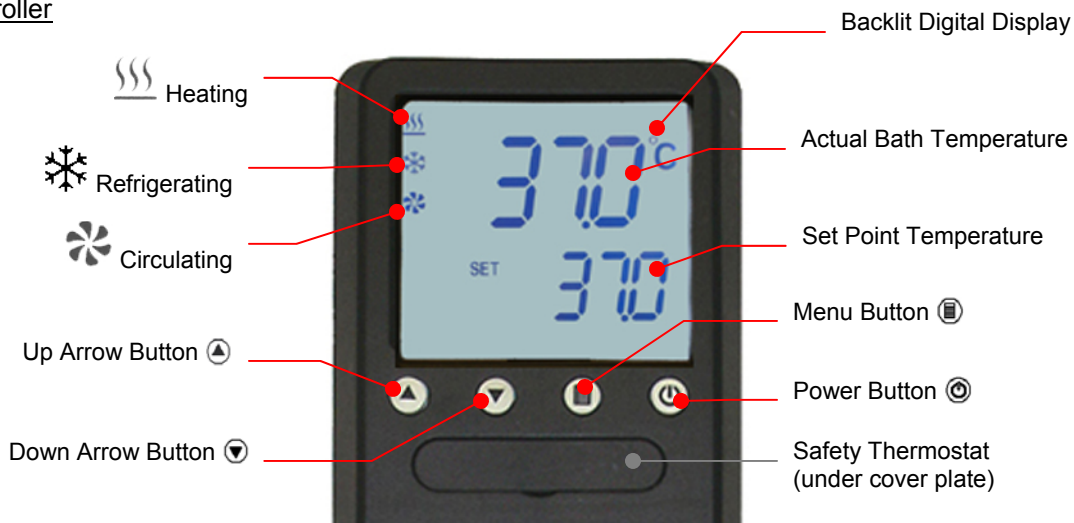

### Refrigerating/Heating Baths



### Open Bath Systems (stainless steel)



**WARNING:** The top deck on Open Bath Systems is not attached. Do not remove deck while Circulator is operating. Do not lift bath by grasping the Temperature Controller or top deck. Always disconnect electrical power and drain fluid from bath before moving.



**WARNING:** To avoid the potential for burns, allow the Circulator to cool completely before cleaning or performing any maintenance.

### Open Bath Systems (polycarbonate)



**CAUTION:** The top deck on Open Bath Systems is not attached. Do not remove deck while Circulator is operating. Do not lift bath by grasping the Temperature Controller or top deck. Always disconnect electrical power and drain fluid from bath before moving.



**WARNING:** To avoid the potential for burns, allow the Circulator to cool completely before cleaning or performing any maintenance.

## Viscosity Baths



**CAUTION:** The top deck on Viscosity Baths is not attached. Do not remove deck while Circulator is operating. Do not lift bath by grasping the Temperature Controller or top deck. Always disconnect electrical power and drain fluid from bath before moving.



**WARNING:** To avoid the potential for burns, allow the Circulator to cool completely before cleaning or performing any maintenance.






## Colform Bath





## Quick-Start

Unless otherwise specified, quick-start instructions apply to all models.

See Installation & Startup for additional information.
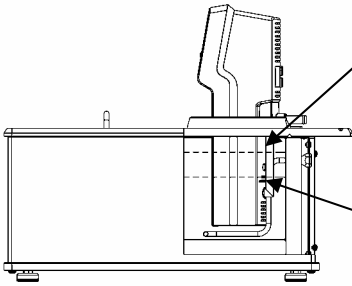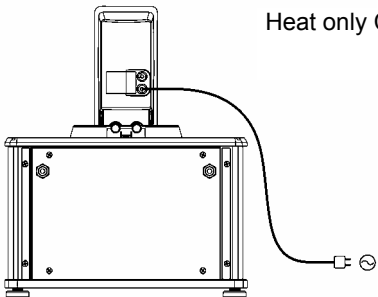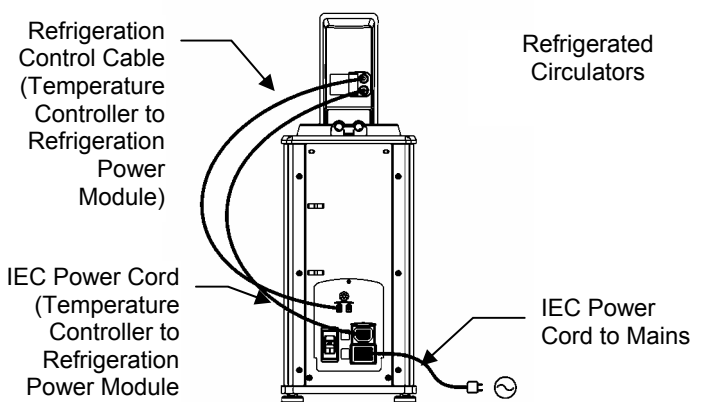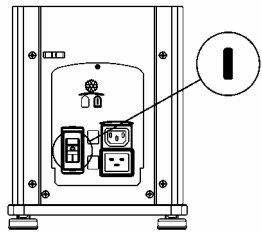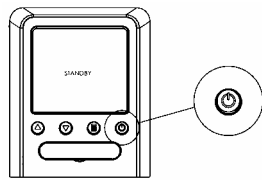
|   |  |   |
|---|--|---|
| 1 | Fill reservoir with fluid  |  <p>Maximum: 1 in / 2.54 cm<br/>below the underside of the top<br/>deck</p> <p>Minimum: 2.6 in. / 6.7 cm<br/>below the underside of the top deck</p>  |
| 2 | Connect all power cords and control cables                                       |  <p>Heat only Circulators</p>   |
|   |  |  <p>Refrigeration Control Cable<br/>(Temperature Controller to Refrigeration<br/>Power Module)</p> <p>IEC Power Cord<br/>(Temperature Controller to Refrigeration<br/>Power Module)</p> <p>Refrigerated Circulators</p> <p>IEC Power Cord to Mains</p> |
| 3 | Turn Power Switch / Circuit Breaker ON<br><br>(Refrigerating/Heating Baths only) |   |
| 4 | Turn Controller ON   |   |
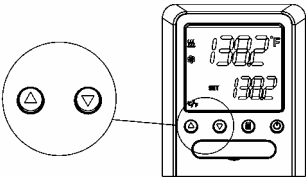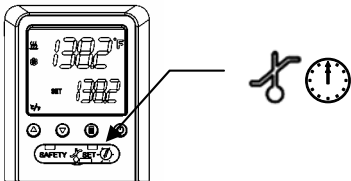
|   |   |  |
|---|---|--|
| 5 | Enter temperature set point                       |  |
| 6 | Set safety thermostat once unit reaches set point |  |

## Installation and Startup

Your Circulating Bath with MX Temperature Controller is designed to be simple to set-up and install. The only tools required are a flat-head screwdriver and a container for adding water or another suitable fluid to the bath reservoir.

### General Site Requirements

Locate your Circulator on a level surface free from drafts and direct sunlight. Do not place it where there are corrosive fumes, excessive moisture, high room temperatures, or in excessively dusty areas.

Refrigerating / Heating Circulators must be 10.2 cm / 4 inches or more away from walls or vertical surfaces so that airflow is not restricted.

Avoid voltage drops by using properly grounded power outlets wired with 14 gauge or larger diameter wire and if possible, be close to the power distribution panel. The use of extension cords is not recommended; this will reduce the potential for problems caused by low line voltage.

### Adding Liquid to the Bath Reservoir




**WARNING:** Read the safety data sheet for the bath fluid being used carefully before filling reservoir.

**WARNING:** Always use fluids that satisfy safety, health, and equipment compatibility requirements.

**WARNING:** If the proper fluid level is not maintained, the heater coil may become exposed and possibly damaged (fluid level too low) or the bath may overflow (fluid level too high).

The liquid in the reservoir should be maintained at a depth between 1 inch / 2.54 cm and 2.6 inches / 6.7 cm below the underside of the bath's top deck. Upon start up, it may be necessary to add fluid to the bath to compensate for the fluid required for external circulation. Likewise, be sure to compensate for fluid displacement when placing samples or other materials in the Circulator's reservoir.

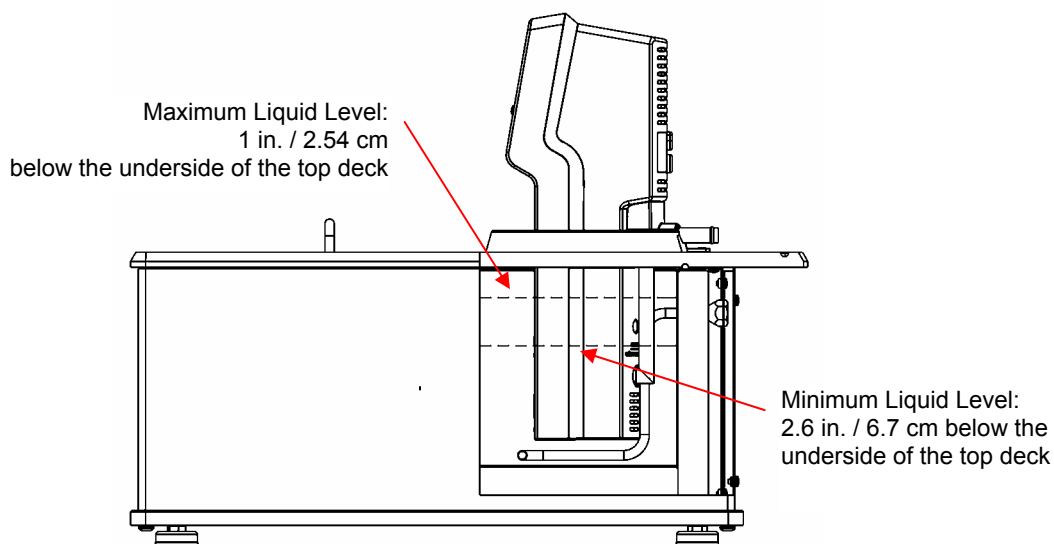



**WARNING:** Always drain all fluid from the reservoir before moving or lifting your Circulator. Be sure to follow your organization's procedures and practices regarding the safe lifting and relocation of heavy objects.



**WARNING:** To avoid the potential for burns, allow the Circulator to cool completely before cleaning or performing any maintenance.

## External Closed Loop Circulation


|   |   |
|---|---|
|  | <p><b>WARNING:</b> When connecting tubing to an external application, it is the user's responsibility to make sure that the tubing and fittings connected to the Circulator are suitable for the fluid being used and the temperature range of operation.</p> <p><b>CAUTION:</b> The Circulator's bypass tubing is secured to the fluid inlet and outlet connections by high temperature nylon hose clamps, which can be removed by carefully cutting them with diagonal cutters.</p> <p><b>CAUTION:</b> Secure the tubing to the inlet and outlet fittings using hose clamps with a minimum ID of 7/8 inch (22 mm). Do not operate the unit without hose clamps.</p> |
|---|---|

|   |   |
|---|---|
|  | <p><b>WARNING:</b> If the Circulating Bath will not be used for external circulation, the inlet and outlet ports should remain connected using the Buna N bypass tubing provided with the unit.</p> |
|---|---|

The single speed pump on your Circulating Bath permits closed loop circulation to an external apparatus via the ½ inch / 12.7 mm OD pump inlet and outlet ports on the rear of the MX Controller. To maintain adequate flow, avoid restrictions in the tubing.


The control stability of a closed loop system is better at the external apparatus than in the immediate vicinity of the heater (provided the apparatus control point represents a constant load and is well insulated). For example, if you circulate through a viscometer at 50°C the temperature variation observed in the reservoir may be ±0.1°C, whereas in the viscometer it may be only ±0.05°C.

Although stability is better at the external apparatus control point, depending on the insulation and length of tubing used, the accuracy of temperature may be slightly different than the temperature indicated in the reservoir.

|  |  |
|--|--|
|  | <p><b>NOTE:</b> Viscosity Baths and the Coliform Bath are not equipped for external circulation.</p> |
|--|--|

## Refrigeration Control Connections (Refrigerating/Heating Circulators only)

Attach the Refrigeration Control Cable connected to the rear of the Temperature Controller to the Refrigeration Control Connection on the Refrigeration Power Module.

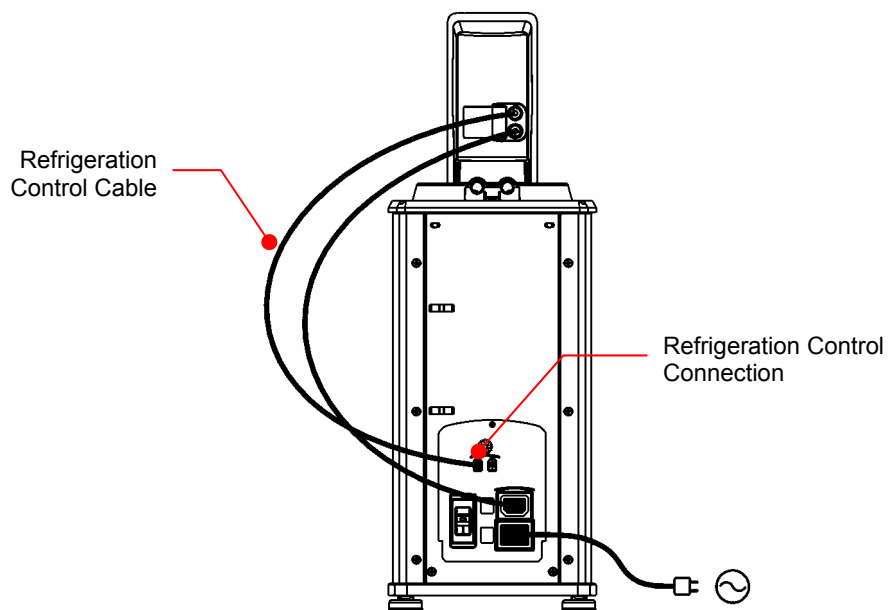



## Electrical Power



**WARNING:** The Circulator's power cord must be connected to a properly grounded electrical receptacle. Make certain that this electrical outlet is the same voltage and frequency as your Circulator. The correct voltage and frequency for your Circulator are indicated on the identification label on the back of the Controller.



**CAUTION:** The use of an extension cord is not recommended. If one is necessary, it must be properly grounded and capable of handling the total wattage of the unit. The extension cord must not cause more than a 10% drop in voltage to the unit.




**NOTE:** All models supplied with European line cords also include line cord extensions for use in the United Kingdom and Switzerland.

### Refrigerating / Heating Circulators

Attach the Temperature Controller's power cord to the female IEC connector on the Refrigeration Power Module.

Attach the 6-ft / 1.8 m power cord to the IEC electrical connection on the Refrigeration Power Module and then plug the male connector into the Mains electrical outlet.

Place the Power Switch / Circuit Breaker on the Refrigeration Power Module in the ON position. The LCD on the Controller will light and "Standby" will appear on the display.

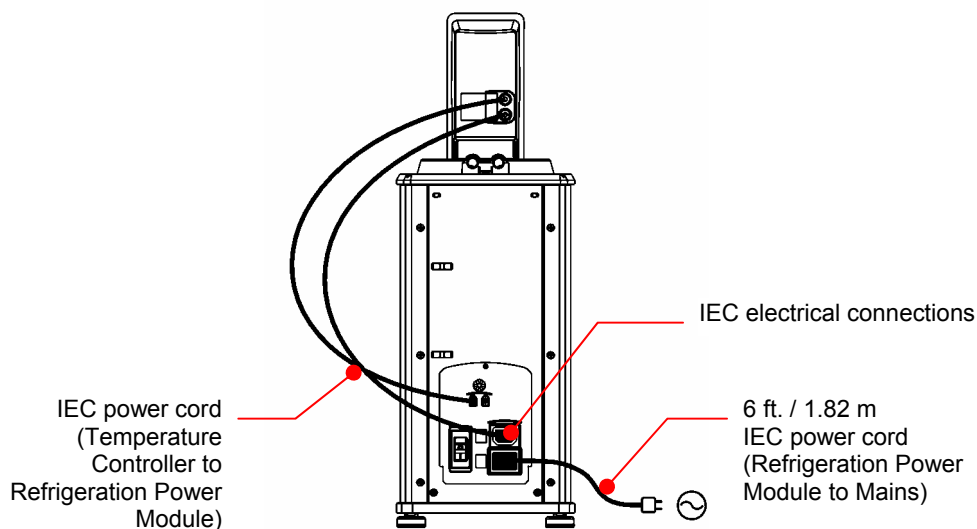


### All Other Circulators

Plug the male connector on the Temperature Controller's power cord into the Mains electrical outlet. The LCD on the Controller will light and "Standby" will appear on the display.




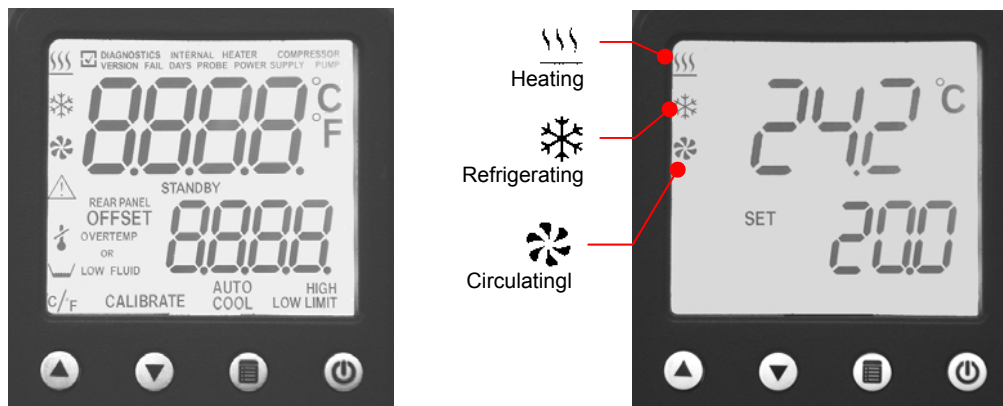
**NOTE:** 50 Hz models supplied with a country specific electrical cord and plug.

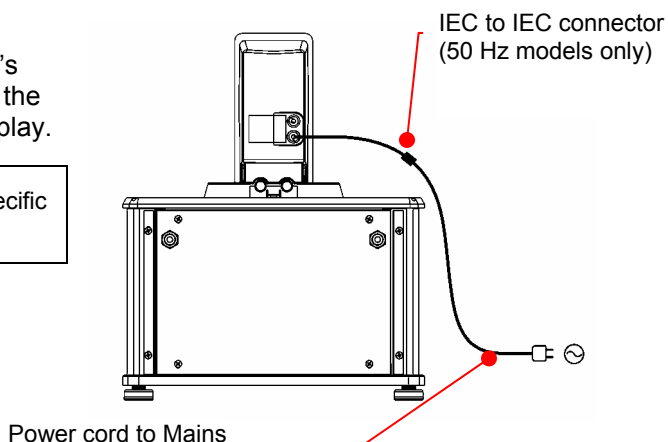

## Controller Setup


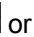


### Power

Press ; all characters and symbols on the LCD will momentarily light. The Circulator will begin running, actual and set point temperatures will be displayed, and the word “SET” will be continuously lit. The circulating symbol will also be lit and the heating or refrigerating symbol may be lit or flashing.



### Safety Set Temperature

This is a “Do Not Exceed” temperature setting for your Circulator and is the temperature at which the heater will be turned OFF should the liquid level in the bath drop too low or the Circulator malfunctions. It is normally set about 5° higher than the desired operating temperature. Setting the Safety Set temperature is a four-step process.

1. Using a flat blade screwdriver, rotate the Safety Thermostat (located beneath the cover plate) clockwise until it stops.
2. Press the  or  key; “SET” will begin flashing. Continue pressing the  and  keys until the set point temperature is equal to your desired Safety Set temperature. “SET” will stop flashing approximately 10 seconds after the desired temperature has been entered. Allow the Circulator to stabilize at this temperature.
3. Once the bath temperature has stabilized, slowly rotate the Safety Thermostat counter-clockwise until you hear a soft “click;” the OVERTEMP or LOW FLUID alarm symbol will appear on the display approximately 5-10 seconds later. At this point, the heater will also turn OFF.
4. Slowly rotate the Safety Thermostat clockwise until the alarm symbol extinguishes (you may also hear a soft “click”). The heater should turn back ON. You are now ready to start normal operation.







OVERTEMP / LOW FLUID Alarm




**WARNING:** The Safety Thermostat is user-adjustable from approximately 40° to 210°C. Do not force the indicator dial beyond the stops at either end of the dial's range.

## Normal Operation

### Keys and Controls

|            |   |  |
|------------|---|--|
| Power      |  | Turns the Circulator's Temperature Controller ON.  |
| Menu       |  | Accesses the Temperature Controller's set-up sub-menus. The items in these sub-menus are used to configure the Controller's general operational parameters (temperature unit, upper and lower temperature limits, offset calibration, etc.). |
| Up Arrow   |  | Used to increase temperature set point and other operational settings/values.  |
| Down Arrow |  | Used to decrease temperature set point and other operational settings/values.  |

### Turning Your Circulator ON

Press the  key.

All characters/symbols on the LCD will momentarily light. When the Circulator begins running, the actual and set point temperatures will be displayed and the circulating symbol will be lit.




If the actual bath temperature is lower than the set point temperature, the heating symbol will also be lit.


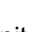

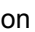






Refrigerating/Heating Models: If the actual bath temperature is higher than the set point temperature, the refrigerating symbol will be lit. It is normal for both the heating and refrigerating symbols to be lit simultaneously when nearing or maintaining the set point temperature.

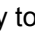
### Main Operational Display



## Set-Up Sub-Menus


Press the  key to access the Temperature Controller's set-up sub-menus. The  or  keys are used to change the current setting / value in the sub-menu.





| Sub-Menu           | Selection / Range               | To Change Current Value  |
|--------------------|---------------------------------|--|
| Temperature unit   | °C or °F                        | Press  for °C<br>Press  for °F |
| Offset Calibration | -3.0°C to +3.0°C                | Press  or                      |
| Low Limit          | -55° to +10°C / -65° to +50°F   | Press  or                      |
| High Limit         | +40° to +155°C / +100 to +305°F | Press  or                      |
| Auto Cool          | 1° to 135°C / 34° to 275°F      | Press  or                      |


To accept a value in a sub-menu, press  or allow the display to 'time out' (approximately 10 seconds).

## Adjusting the Temperature Set Point

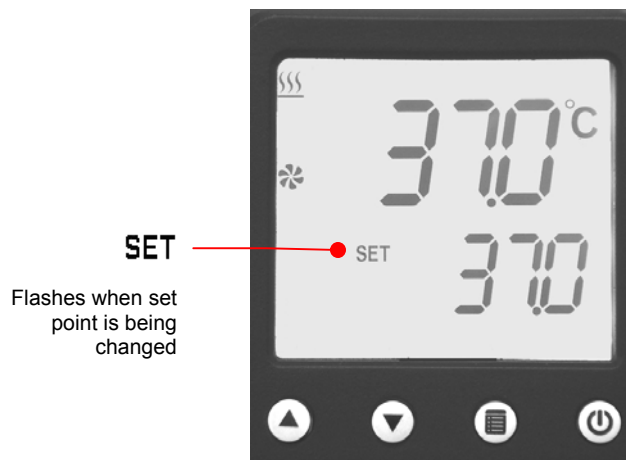This is the temperature at which the fluid in your Circulating Bath will be maintained. It may be set to one-tenth of a degree over a range of -50.0° to +135.0°C / -60.0° to +275.0°F. The factory default set point is +20.0°C / +68.0°F.

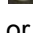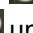



**To Change:** Press the  or  key. The word "SET" will begin flashing. Continue holding or repeatedly press  or  until the desired set point temperature is displayed.

**To Accept:** Press  or allow the LCD to return to the main operational display (approximately 10 seconds).


## Selecting the Temperature Unit

The temperature units sub-menu (°C / °F) allows you to select the temperature unit in which the actual bath temperature and set point temperature are displayed. The factory default is °C.

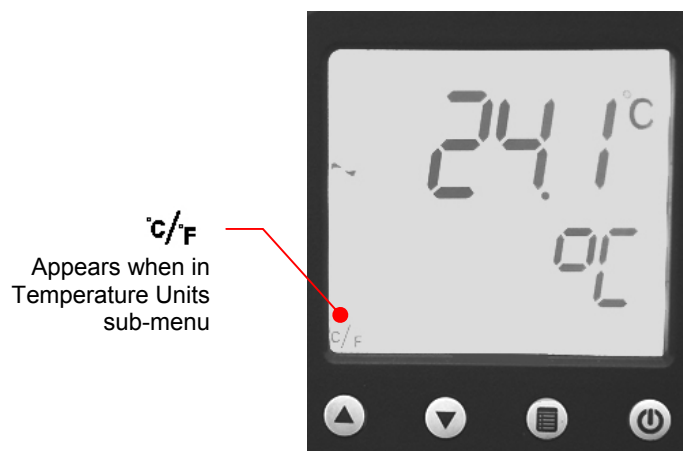








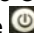
**To Access:** Press the  key until °C/°F appears on the display.


**To Change:** Press  to select °C; press  to select °F.

**To Accept:** Press  or allow the LCD to return to the main operational display (approximately 10 seconds).

## Setting the Offset Calibration


This sub-menu allows you to match the Circulator's temperature display to an external reference thermometer. A value from -3.0° to +3.0°C may be entered; the factory default is 0.0°C.


|   |   |
|---|---|
|  | <b>IMPORTANT:</b> To prevent the Offset Calibration value from being changed unintentionally, the following power down/power up sequence is required to enable the Offset Calibration function.   |
|   | <p><u>Refrigerating Circulators:</u></p> <ol style="list-style-type: none"><li>1. Place the power switch/circuit breaker on the rear of the unit in the OFF position.</li><li>2. Return the power switch/circuit breaker to the ON position while pressing and holding the  key.</li><li>3. When STANDBY appears on the display, release the  key and press . You can now proceed as outlined below.</li></ol> <p><u>Heat Only Circulators:</u></p> <ol style="list-style-type: none"><li>1. Disconnect the power cord from the electrical outlet.</li><li>2. Plug the power cord back into the electrical outlet while pressing and holding the  key.</li><li>3. When STANDBY appears on the display, release the  key and press . You can now proceed as outlined below.</li></ol> <p>The Offset Calibration function will remain enabled until the electrical power is turned OFF by pressing the  key.</p> |

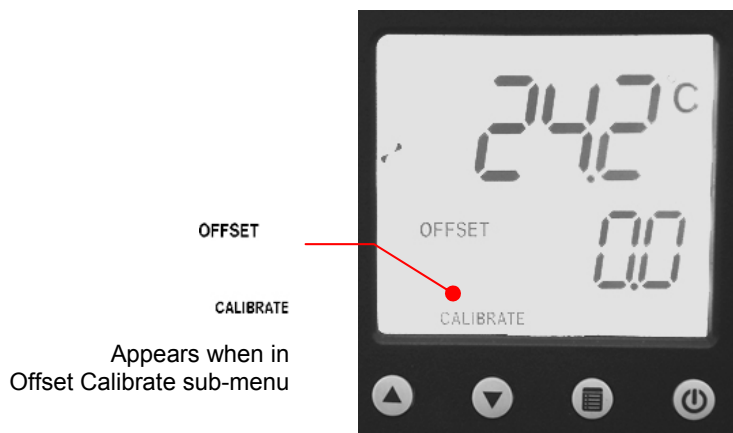

**To Access:** Press the  key until OFFSET CALIBRATE appears on the display.

**To Change:** Press  or .

**To Accept:** Press  or allow the LCD to return to the main operational display (approximately 10 seconds).

|   |  |
|---|--|
|  | <b>CAUTION:</b> The Offset Calibration value is always shown in degrees C, even if degrees F is selected as the temperature unit in which the control and actual bath temperatures are displayed. Your Circulator will automatically convert the °C offset calibration value to the correct °F display offset value. |
|---|--|

## Setting the Low Limit Temperature

This sub-menu allows you to limit how low the temperature set point may be set. It also serves as a low limit safety, alerting you if bath temperature falls below the low limit temperature setting. The Low Limit value may be set from -55° to +10°C / -65° to +50°F; the factory default is -50°C / -60°F.

To avoid unwanted alarms or shutdowns during regular operation, the Low Limit value should be set at least 5° lower than the selected control temperature.




**LOW LIMIT**

Appears when in the Low Limit sub-menu

**To Access:** Press the  key until LOW LIMIT appears on the display.

**To Change:** Press  or .

**To Accept:** Press  or allow the LCD to return to the main operational display (approximately 10 seconds).

## Setting the High Limit Temperature

This sub-menu allows you to limit how high the temperature set point may be set. It also serves as a high limit safety, alerting you if bath temperature rises above the high limit temperature setting. The High Limit value may be set from +40° to +155°C / +100° to +305°F; the factory default is +150°C / +300°F.

To avoid an unwanted shutdown during regular operation, the High Limit value should be set at least 5° higher than the selected control temperature.




**HIGH LIMIT**

Appears when in the High Limit sub-menu

**To Access:** Press the  key until HIGH LIMIT appears on the display.

**To Change:** Press  or .

**To Accept:** Press  or allow the LCD to return to the main operational display (approximately 10 seconds).

## Adjusting the Auto Cool Temperature

This sub-menu is displayed only on Refrigerating / Heating Circulators. It determines the bath temperature at which refrigeration will be activated and permits more precise control when operating at high temperatures as well as more rapid cool downs. For most applications, an Auto Cool set point 15°C above room temperature is recommended. The Auto Cool range is from 1° to 135°C. The factory default is 30°C.


Cool Command™ Refrigeration — Refrigerating/Heating Circulators with 15 liter and larger reservoirs feature the Cool Command™ modulating refrigeration control system. Cool Command™ allows the refrigeration system to turn on at a fluid temperature up to 135°C when the temperature set point is changed to or below the Auto Cool set point (135°C maximum). As a result, bath fluid cools more quickly.

Conventional Refrigeration — Refrigerating/Heating Circulators with 7 liter reservoirs use a conventional refrigeration system. The refrigeration system will turn on when the bath temperature and set point are below the Auto Cool set point (85°C maximum).




**AUTO  
COOL**

Appears when in the  
AUTO COOL menu

**To Access:** This Press the  key until AUTO COOL appears on the display.




**To Change:** Press  or .

**To Accept:** Press  or allow the LCD to return to the main operational display (approximately 10 seconds).



## Resetting the Factory Default Values

You can return your Circulator to all factory-set defaults as follows:

1. Press the  key to turn the Circulator OFF.
2. Turn Mains power OFF by unplugging the power cord from the electrical outlet or (Refrigerating/Heating Circulators) placing the power switch/circuit breaker on the Refrigeration Power Module in the OFF position.
3. Turn Mains power ON while pressing either the  or  button.


The factory default values are:

| Operational Parameter | Factory Default Value |
|-----------------------|-----------------------|
| Temperature Scale     | °C                    |
| Temperature Set Point | 20°C / 68°F           |
| Calibration Offset    | 0.0°C                 |
| Low Limit             | -50°C / -60°F         |
| High Limit            | 150°C / 300°F         |
| Auto Cool             | 30°C                  |

## Loss of Power Restart



**WARNING:** The unit will start automatically after a disruption in electrical power.

In the event that electrical power is lost while your Circulator is in use, it will begin operating automatically once power is restored. FAIL POWER will appear on the display to alert you that there was a power disruption. To clear the message, turn the Circulator OFF and then back ON again using the  key.



**FAIL POWER**

Indicates electrical power was disrupted during operation

## Tap Water Cooling

Tap water cooling allows for more rapid bath cool down from high temperatures and/or more precise operation at temperatures near ambient.

Viscosity Baths feature a cooling coil with two 0.375 inch / 9.5 mm OD straight barb fittings at which the fluid inlet and outlet connections can be made. Either connection may serve as the fluid inlet or outlet. Be sure to secure the tubing with the appropriate size hose clamps.


An optional cooling coil is available for use on the Coliform Bath, Stainless Steel Open Bath Systems, and Polycarbonate Open Bath Systems. This cooling coil has two 0.375 inch / 9.5 mm OD straight barb fittings at which the fluid inlet and outlet connections can be made. Either connection may serve as the fluid inlet or outlet. Be sure to secure the tubing with the appropriate size hose clamps.







**WARNING:** The fluid outlet must be connected and flow to a suitable drain or vessel located at a level below that of the inlet.
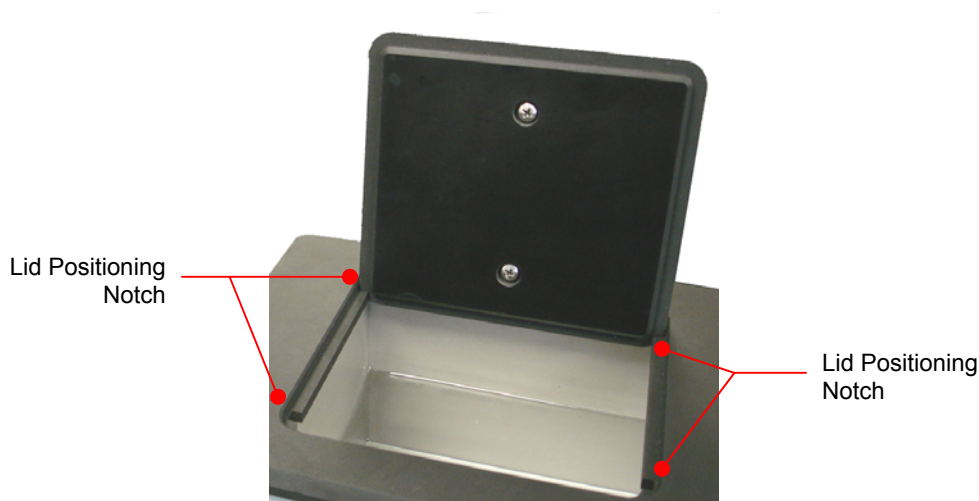
## Reservoir Cover Storage

Refrigerating/Heating and Heat Only Circulating Baths feature the LidDock<sup>®</sup> system to eliminate mess when adding fluid or samples to the reservoir. Specially positioned notches in the inner lip of the top deck allow you to stand the reservoir cover up upright over the bath opening, allowing condensate to flow back into the bath.



## Display Messages and Alarms

| Alarm Symbol  | Description  | Corrective Action  |
|---|--|--|
| <b>FAIL POWER</b>   | <b>Informational Message:</b><br>Indicates that electrical power was lost during operation.  | Using the  key, turn the Circulator OFF and then back ON. This will clear the message.  |
| <b>LOW LIMIT</b><br>(flashing)  | <b>Warning:</b> The temperature set point is below the Low Limit temperature value.  | Decrease the Low Limit temperature value or increase the temperature set point.  |
| <b>HIGH LIMIT</b><br>(flashing)   | <b>Warning:</b> The temperature set point is above the High Limit temperature value.   | Increase the High Limit temperature value or decrease the set point temperature.   |
| <br><b>LOW LIMIT</b>               | <b>Alarm:</b> The bath temperature has fallen below the Low Limit temperature value. Power to the compressor and pump will remain OFF until the problem is corrected.                                      | Allow bath to warm or add heat load.<br>Decrease the Low Limit temperature value.  |
| <br><b>HIGH LIMIT</b>              | <b>Alarm:</b> The bath temperature has risen above the High Limit temperature value. Power to the heater and pump will remain OFF until the problem is corrected.  | Allow bath to cool or increase High Limit temperature value.<br>Replace fluid.   |
| <br><b>OVERTEMP OR LOW FLUID</b> | <b>Fault:</b> The liquid in the bath has dropped too low or the temperature of the bath fluid has exceeded the Safety Set temperature. Power to the heater will remain OFF until the problem is corrected. | Fluid level in reservoir has fallen below minimum level; add fluid as required.<br>Fluid temperature is higher than Safety Set temperature; increase Safety Set temperature setting.<br>Controller failure; consult factory. |
| <b>INTERNAL PROBE FAIL</b>  | <b>Fault:</b> The Circulator's temperature sensor has failed.  | Consult factory.   |
| <b>HEATER FAIL</b>  | <b>Fault:</b> The Circulator's heater has failed.  | Consult factory.   |

## Routine Maintenance and Troubleshooting



**WARNING:** Always turn your Circulator OFF and disconnect it from the electrical power outlet before performing any maintenance or service.



**WARNING:** To avoid the potential for burns, allow the Circulating Bath to cool completely before cleaning or performing any maintenance.



**WARNING:** Always drain all fluid from the reservoir before moving or lifting your Circulator. Be sure to follow your organization's procedures and practices regarding the safe lifting and relocation of heavy objects.

### Maintaining Clear Bath Water

Optimum temperature and moisture conditions for algae growth exist when using water as a bath fluid. To prevent algae contamination and minimize the frequency of draining the reservoir, an algaecide should be used.



**WARNING:** Do not use chlorine bleach.

## Draining the Bath Reservoir




**WARNING:** Bath fluids should be stored and disposed of according to applicable laws and regulations.

### Refrigerated / Heating Circulating Baths

Refrigerated / Heating and Heat only Circulating Baths are equipped with a drain valve and port located either behind the front access panel or on the right hand side of the unit.

To drain fluid from the bath, attach a short length of suitable 11.5 mm ID / 0.45 inch ID tubing to the drain port and secure it using a hose clamp with a minimum ID of 18 mm / 0.7 inch. Open the drain valve using a flat blade screwdriver. When closing the valve, do not over tighten.

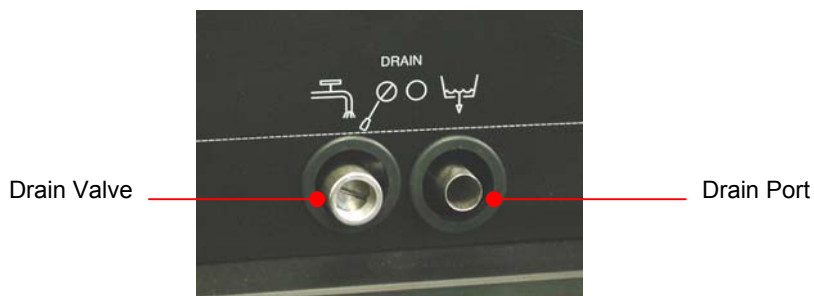

**WARNING:** Be sure to close the drain valve before refilling the bath reservoir. Do not over tighten.

### Polycarbonate Open Tank Circulators and Viscosity Baths

Larger capacity (17 liter and above) polycarbonate open tank Circulating Baths and Viscosity Baths are equipped with a drain port. To drain fluid from these baths, remove the knurled cap. When replacing the cap, do not over tighten.



**WARNING:** Be sure to replace and tighten knurled cap before refilling the bath reservoir

## Checking the Over-Temperature Safety System

Your Circulator incorporates over-temperature protection according to IEC 61010. For optimum safety, this system should be checked for proper operation at least every six months. This check must be performed with the unit running.

1. Enter a temperature set point of approximately 50°C and then allow the bath to stabilize at that temperature. The amount of time this will take will depend on the size of the bath and the difference between the initial bath temperature and the Safety Set temperature.
2. Once the bath temperature has stabilized, slowly rotate the Safety Thermostat counter-clockwise until you hear a soft “click;” the OVERTEMP or LOW FLUID alarm symbol will appear on the display approximately 5-10 seconds later. At this point, the heater will also turn OFF.
3. Slowly rotate the Safety Thermostat clockwise until the alarm symbol extinguishes (you may also hear a soft “click”). The heater should turn back ON.
4. Reset the Safety Set temperature to the desired temperature value (see *Controller Setup, Safety Set Temperature*).



## Cleaning Your Circulator



**WARNING:** It is the user's responsibility to properly decontaminate the unit in the event hazardous materials are spilled on exterior or interior surfaces. Consult the manufacturer if there is any doubt regarding the compatibility of decontamination or cleaning agents.

### Temperature Controller

Turn the Temperature Controller OFF by pressing  and unplug power cord from the electrical outlet.

Wipe the housing with a clean cloth dampened with a mild detergent and water or mild all-purpose cleaner.



**CAUTION:** Do not spray cleaning liquids directly onto the Temperature Controller or allow them to enter the Controller's vents. Do not use abrasives as these could scratch the housing or the digital display.

### Bath Reservoir



**CAUTION:** Do not use steel wool to clean your Circulator's bath reservoir.

**Bath Reservoir and Wetted Components** — A vinegar solution or calcium/lime/rust remover can be used to clean mineral deposits from the Temperature Controller's wetted parts. The cleaner should be added to the bath reservoir at the prescribed dosage and circulated at 60°C / 140°F until the scale is removed.

**External Surfaces** — Only mild detergents and water or an approved cleaner should be used on the top deck and other external surfaces of your Circulator.

**Pump Impeller** — In the unlikely event that debris becomes lodged in the pump impeller, a soft brush can be used to remove any lodged particles. If necessary, soak in a solution of distilled water and vinegar to soften before brushing.




**CAUTION:** Do not use hard utensils or abrasive pads to remove trapped debris.

### Condenser, Air Vents, and Reusable Filter (Refrigerated Models only)

To keep the refrigeration system operating at optimum cooling capacity, the condenser, removable air filter, and all air vents (front, side, back) should be kept free of dust and dirt. Be sure to check them on a regular basis and clean as required.

The reusable filter is easily accessed from the front of the unit by simply removing the access panel. Use a mild detergent and water solution to wash off any accumulated dust and dirt. Rinse and dry thoroughly before reinstalling.

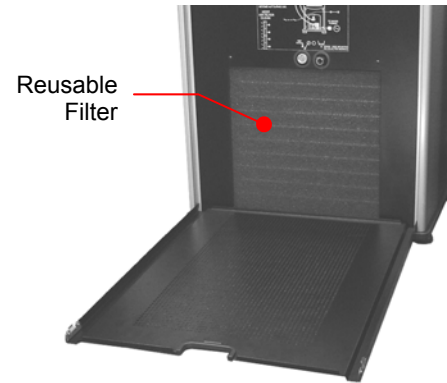


## Temperature Controller Removal and Re-Installation


### Removal

The Temperature Controller on your Circulating Bath is designed to be easily removed from the top deck without the use of special tools. It is removed as follows:


1. Place the tip of a small flat blade screwdriver under the retaining ring locking tab and pry up gently.

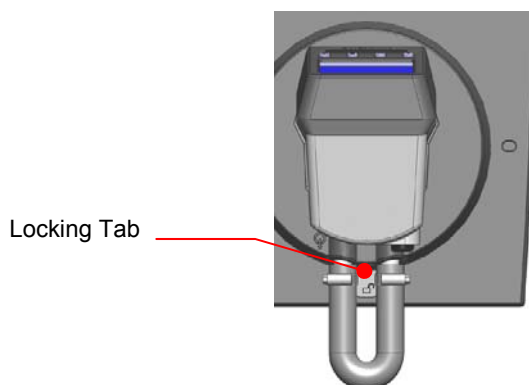

2. Rotate the Temperature Controller clockwise until it stops (about 0.75 inch / 1.9 cm).

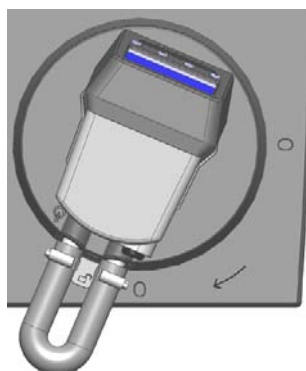

3. Lift the Controller straight up and out of the opening of the Circulator's top deck.

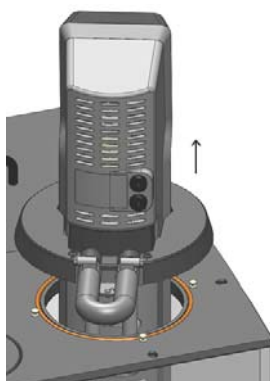





## Re-Installation

The top deck of your Circulator incorporates four alignment pins to facilitate positioning of the Temperature Controller when it is being reinstalled. These pins correspond to keyhole slots on the interior of the Circulator's retaining ring.


1. With the retaining ring locking tab oriented above one of the indents on the top deck, slowly lower the Temperature Controller into the top deck opening until it is resting on top of the positioning pins.

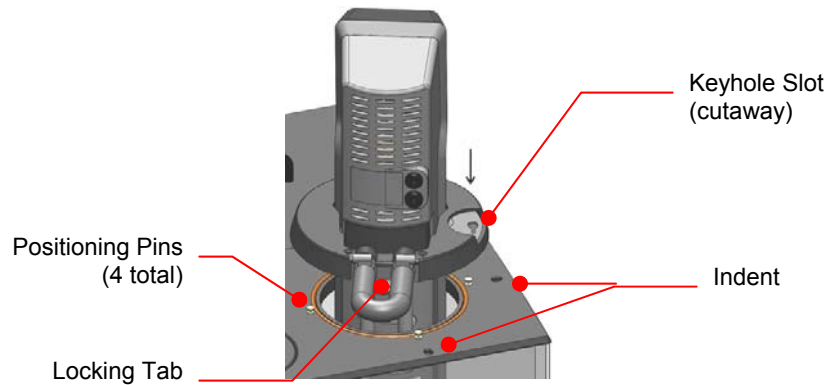

2. Gently rotate the Temperature Controller until it drops down on the Positioning Pins.

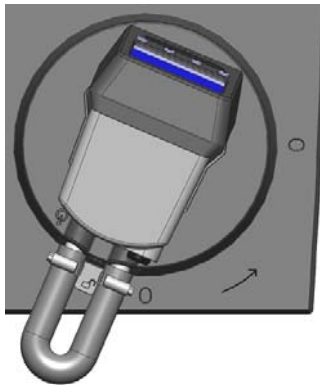

3. Rotate the Temperature Controller counterclockwise until the Locking Tab engages the indent on the top deck.

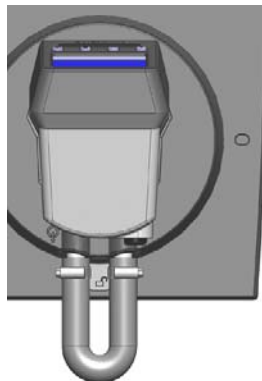

## Troubleshooting Chart

| Problem   | Possible Causes  | Corrective Action  |
|---|--|--|
| Unit does not run<br>(Digital Display is blank)           | No power to unit<br><br><i>Refrigerated units only</i><br>Electrical cord between Control Head and Refrigeration Power Module loose or unplugged<br>Power Switch / Circuit Breaker in OFF position                         | Check that the electrical cord is secure and connected to an operating electrical outlet.<br><br>Check that the electrical cord is secure and properly connected.<br><br>Place Power Switch / Circuit Breaker in ON position.  |
| Unit does not run<br>(STANDBY appears on Digital Display) | Unit in Standby mode   | Press Power Key on front panel.  |
| No fluid circulation                                      | Insufficient fluid in reservoir<br>Pump impeller jammed  | Add fluid to reservoir.<br>Inspect pump and remove debris as required.   |
| Insufficient circulation                                  | Fluid viscosity too high<br>External tubing diameter too small<br>Low line voltage   | Replace with lower viscosity bath fluid.<br>Replace with larger diameter tubing.<br><br>Check and correct as required.   |
| Unit does not heat  | Insufficient fluid in reservoir<br>Temperature set point too low<br>Safety Set Temperature too low   | Add fluid to reservoir.<br>Increase temperature set point.<br><br>Increase Safety Set temperature.   |
| Insufficient heating                                      | Insufficient circulation<br>Low line voltage<br>Ambient temperature too cool<br>Excessive heat loss  | See Insufficient circulation, above.<br>Check and correct as required.<br>Increase ambient temperature or relocate unit.<br><br>Check for heat loss from external tanks and hoses;<br>Check for vapor/heat loss from internal reservoir.   |
| Temperature unstable                                      | Insufficient circulation<br>Debris or mineral build-up on pump, heater, or temperature sensor.   | Check pump flow and operation.<br>Clean as required.   |
| Unit does not cool  | Dust build up on air filter or condenser<br>Blocked air ventilation screens<br>Temperature set point is too high<br>Excessive heat load<br><br>Ambient air temperature too high (>35°C / 95°F)<br>Low or high line voltage | Clean air filter and/or condenser as required.<br><br>Remove blockages as required.<br><br>Decrease temperature set point.<br><br>Check that heat load does not exceed capacity of bath; correct as required.<br>Decrease ambient air temperature.<br><br>Check and correct as required. |

| <b>Problem</b>                                 | <b>Possible Causes</b>  | <b>Corrective Action</b>   |
|--|---|--|
| Insufficient cooling                           | Dust build up on air filter or condenser<br>Blocked air ventilation screens<br>Temperature set point is too high<br>Excessive heat load<br>Ambient air temperature too high (>35°C / 95°F)<br>Low or high line voltage  | Clean air filter and/or condenser as required.<br>Remove blockages as required.<br>Decrease temperature set point.<br>Check that heat load does not exceed capacity of bath; correct as required.<br>Decrease ambient air temperature.<br>Check and correct as required.   |
| Unable to achieve low end extreme temperatures | Pump speed too high<br>Incorrect bath fluid<br>Insufficient insulation on external fluid lines<br>Ambient air temperature too high (>35°C / 95°F)<br>Low or high line voltage<br>Dust build up on air filter or condenser<br>Blocked air ventilation screens<br>Excessive heat load | Reduce pump speed.<br>Check that the fluid being circulated is capable of reaching the required temperature.<br>Check external fluid lines for proper insulation.<br>Decrease ambient air temperature as required.<br>Check and correct as required.<br>Clean air filter or condenser as required.<br>Remove blockages as required.<br>Check that heat load does not exceed capacity of bath; correct as required. |

# Technical Information

## Performance Specifications

|                              |                                  |                    |
|------------------------------|----------------------------------|--------------------|
| Operating Temperature Range: | Model dependent; see table below |                    |
| Temperature Stability:       | ±0.07C (±0.13°F)                 |                    |
| Pump Type:                   | 1-speed pressure                 |                    |
|                              | <u>60Hz models</u>               | <u>50Hz models</u> |
| Maximum Pressure:            | 1.8 psi (0.12 bar)               | 1.5 psi (0.10 bar) |
| Maximum Pressure Flow Rate:  | 3.4 gpm (12.8 lpm)               | 2.8 gpm (10.6 lpm) |
| Heater Wattage:              | 1100 watts                       | 1100 watts         |

| Model Type                              | Reservoir Capacity | Temperature Range  | Electrical Requirements |                |
|---|--------------------|--|-------------------------|----------------|
|   |                    |  | 60Hz Units              | 50Hz Units     |
| MX07R-20 Refrigerating / Heating Bath   | 7 liters           | -20° to 135°C<br>-4° to 275°F                                | 120V, 60Hz, 12A         | 240V, 50Hz, 8A |
| MX7LR-20 Refrigerating / Heating Bath   | 7 liters           | -20° to 135°C<br>-4° to 275°F                                | 120V, 60Hz, 12A         | 240V, 50Hz, 8A |
| MX06S135 Stainless Steel Open Tank Bath | 6 liters           | Ambient +10° to 135°C<br>Ambient +20° to 275°F               | 120V, 60Hz, 10A         | 240V, 50Hz, 6A |
| MX10S135 Stainless Steel Open Tank Bath | 10 liters          | Ambient +10° to 135°C<br>Ambient +20° to 275°F               | 120V, 60Hz, 10A         | 240V, 50Hz, 6A |
| MX20S135 Stainless Steel Open Tank Bath | 20 liters          | Ambient +10° to 135°C<br>Ambient +20° to 275°F               | 120V, 60Hz, 10A         | 240V, 50Hz, 6A |
| MX11P100 Polycarbonate Open Tank Bath   | 11 liters          | Ambient +10° to 85°C<br>Ambient +20° to 185°F <sup>(1)</sup> | 120V, 60Hz, 10A         | 240V, 50Hz, 6A |
| MX17P100 Polycarbonate Open Tank Bath   | 17 liters          | Ambient +10° to 85°C<br>Ambient +20° to 185°F <sup>(1)</sup> | 120V, 60Hz, 10A         | 240V, 50Hz, 6A |
| MX28P100 Polycarbonate Open Tank Bath   | 28 liters          | Ambient +10° to 85°C<br>Ambient +20° to 185°F <sup>(1)</sup> | 120V, 60Hz, 10A         | 240V, 50Hz, 6A |
| MX29VB3S Viscosity Bath                 | 29 liters          | Ambient +10° to 85°C<br>Ambient +20° to 185°F <sup>(1)</sup> | 120V, 60Hz, 10A         | 240V, 50Hz, 6A |
| MX29VB5R Viscosity Bath                 | 29 liters          | Ambient +10° to 85°C<br>Ambient +20° to 185°F <sup>(1)</sup> | 120V, 60Hz, 10A         | 240V, 50Hz, 6A |
| MX28C135 Coliform Bath                  | 28 liters          | Ambient +10° to 135°C<br>Ambient +20° to 275°F               | 120V, 60Hz, 10A         | 240V, 50Hz, 6A |

1. Maximum operating temperature for polycarbonate tank; MX Controller is capable of higher temperatures.

|                          |   |
|--------------------------|---|
| Environmental Conditions | Indoor use only                             |
|                          | Maximum Altitude: 2000 meter                |
|                          | Operating Ambient: 5° to 35°C (41° to 95°F) |
|                          | Relative Humidity: 80%, non-condensing      |
|                          | Installation Category: II                   |
|                          | Pollution Degree: 2                         |
|                          | Ingress Protection: IP 31                   |
|                          | Climate Class: SN                           |
|                          | Software Class: B                           |
|                          | Output Waveform: Sinusoidal                 |

Specifications subject to change without notice.

## Reservoir Fluids

Depending on your needs, a variety of fluids can be used with your Circulator. No matter what bath fluid is selected, it must be chemically compatible with the reservoir and the materials in your Circulator. It must also be suitable for the desired temperature range.



**WARNING:** Do not use a flammable liquid as a bath fluid as a fire hazard may result.



**WARNING:** Always use fluids that satisfy safety, health, and equipment compatibility requirements.

For optimum temperature stability, the fluid's viscosity should be 50 centistokes (cSt) or less at its lowest operating temperature. This permits good fluid circulation and minimizes heating from the pump.

For temperatures from 10°C to 90°C, distilled water is recommended. For temperatures below 10°C, a mixture of laboratory grade ethylene glycol and water should be used. Do not use deionized water.



**WARNING: DO NOT USE THE FOLLOWING LIQUIDS**

- Automotive antifreeze with additives\*\*
- Hard tap water\*\*
- Deionized water with a specific resistance > 1 meg ohm
- Any flammable fluids
- Concentrations of acids or bases
- Solutions with halides: chlorides, fluorides, bromides, iodides or sulfur
- Bleach (Sodium Hypochlorite)
- Solutions with chromates or chromium salts
- Glycerine
- Syltherm fluids

\*\* At temperatures above 40°C, additives or mineral deposits can adhere to the heater. If deposits are allowed to build up, the heater may overheat and fail. Higher temperatures and higher concentrations of additives will hasten deposit build up.

## Application Notes

At a fluid's low temperature extreme:

- The presence of ice or slush adversely affects temperature stability.
- A viscosity above 10 centistokes adversely affects temperature uniformity.
- A high fluid viscosity and high pump speed adds heat to the fluid being pumped.

At a fluid's temperature above ambient without refrigeration:

- If your set point temperature is less than 15°C above the ambient temperature, the viscosity of the fluid should be 10 centistokes or less to minimize friction heating of the fluid.
- Heat loss should be encouraged by uncovering the fluid and lowering the pump speed.

At fluid's high temperature extreme:

- Heat loss from vapor adversely affects temperature stability.
- To prevent the accumulation of vapors inside the room, the reservoir may need to be placed in a fume hood.
- Use a cover and/or floating hollow balls to help prevent heat and vapor loss.
- Replenish fluid lost from vapor frequently.

## Equipment Disposal (WEEE Directive)



or



This equipment is marked with the crossed out wheeled bin symbol to indicate it is covered by the Waste Electrical and Electronic Equipment (WEEE) Directive and is not to be disposed of as unsorted municipal waste. **Any products marked with this symbol must be collected separately, according to the regulatory guidelines in your area.**

It is your responsibility to correctly dispose of this equipment at lifecycle-end by handing it over to an authorized facility for separate collection and recycling. It is also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect the persons involved in the disposal and recycling of the equipment from health hazards. By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Requirements for waste collection, reuse, recycling, and recovery programs vary by regulatory authority at your location. Contact your local responsible body (e.g., your laboratory manager) or authorized representative for information regarding applicable disposal regulations.

## Technical Service

### Web Resources

Visit the VWR's website at [www.vwr.com](http://www.vwr.com) for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

**Contact us** For information or technical assistance contact your local VWR representative or visit. [www.vwr.com](http://www.vwr.com).

## Warranty

**VWR International** warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of purchase. If a defect is present, VWR will, at its option, repair, replace, or refund the purchase price of this product at no charge to you, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear.

For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

## Local VWR Offices in Europe and Asia Pacific

### Austria

VWR International GmbH  
Graumannsgasse 7  
1150 Wien  
Tel.: 01 97 002 0  
Fax: 01 97 002 600  
E-mail: info@at.vwr.com

### Belgium

VWR International bvba  
Researchpark Haasrode 2020  
Geldenaaksebaan 464  
3001 Leuven  
Tel.: 016 385 011  
Fax: 016 385 385  
E-mail:  
customerservice@be.vwr.com

### China

VWR International China Co., Ltd  
Suite 1802 - 1803,  
Xing Ye Bank Mansion, No 168,  
168 Jiangning Road  
Shanghai 200041, China  
Tel.: +86- 21 521 388 22  
Fax: +86- 21 521 33 933  
E-mail: sales\_china@vwr.com

### Denmark

VWR - Bie & Berntsen  
Transformervej 8  
2730 Herlev  
Tel.: 43 86 87 88  
Fax: 43 86 87 90  
E-mail: info@dk.vwr.com

### Finland

VWR International Oy  
Valimotie 9  
00380 Helsinki  
Tel.: +358 9 80 45 51  
Fax: +358 9 80 45 52 00  
E-mail: info@fi.vwr.com

### France

VWR International S.A.S.  
Le Périgares – Bâtiment B  
201, rue Carnot  
94126 Fontenay-sous-Bois cedex  
Tel.: 0 825 02 30 30 (0,15 EUR  
TTC/min)  
Fax: 0 825 02 30 35 (0,15 EUR  
TTC/min)  
E-mail: info@fr.vwr.com

### Germany

VWR International GmbH  
Hilpertstrasse 20a  
D - 64295 Darmstadt  
Tel.: 0180 570 20 00\*  
Fax: 0180 570 22 22\*  
E-mail: info@de.vwr.com  
\*0,14 €/Min. aus d. dt. Festnetz,  
Mobilfunk max. 0,42 €/Min.

### Hungary

VWR Spektrum-3D Kft.  
Simon László u. 4.  
4034 Debrecen  
Tel.: (52) 521-131  
Fax: (52) 470-069  
E-mail: info@spektrum-3d.hu

### India

VWR Lab Products Pte Ltd  
2nd Floor, Front Wing, 135/12,  
Brigade Towers  
Brigade Road  
Bangaluru 560025 India  
Tel: +91-2522-277876 (Mumbai)  
Tel: +91-80-41117124 (Bangalore)  
Fax +91-80-41117120  
E-mail: vwr\_india@vwr.com

### Ireland / Northern Ireland

VWR International Ltd / VWR  
International (Northern Ireland) Ltd  
Orion Business Campus  
Northwest Business Park  
Ballycoolin  
Dublin 15  
Tel.: 01 88 22 222  
Fax: 01 88 22 333  
E-mail sales@ie.vwr.com

### Italy

VWR International s.r.l.  
Via Stephenson 94  
20157 Milano (MI)  
Tel.: 02 332 03 11  
Fax: 800 152 999  
E-mail: info@it.vwr.com

### The Netherlands

VWR International B.V.  
Postbus 8198  
1005 AD Amsterdam  
Tel.: 020 4808 400  
Fax: 020 4808 480  
E-mail: info@nl.vwr.com

### Norway

VWR International AS  
Haavard Martinsens vei 30  
0978 Oslo  
Tel.: 02290  
Fax: 815 00 940  
E-mail: info@no.vwr.com

### Poland

Labart Sp. z o.o.  
A VWR International Company  
Limbowa 5  
80-175 Gdansk  
Tel.: 058 32 38 210  
Fax: 058 32 38 205  
E-mail: labart@labart.pl

### Portugal

VWR International - Material de  
Laboratório, Lda  
Edifício Neopark  
Av. Tomás Ribeiro, 43- 3 D  
2790-221 Carnaxide  
Tel.: 21 3600 770  
Fax: 21 3600 798/9  
E-mail: info@pt.vwr.com

### Singapore

VWR Singapore Pte Ltd  
18 Gul Drive  
Singapore 629468  
Tel: +65 6505 0760  
Fax: +65 6264 3780  
E-mail: sales@sg.vwr.com

### Spain

VWR International Eurolab S.L.  
C/ Tecnología 5-17  
A-7 Llinars Park  
08450 - Llinars del Vallès  
Barcelona  
Tel.: 902 222 897  
Fax: 902 430 657  
E-mail: info@es.vwr.com

### Sweden

VWR International AB  
Fagerstagatan 18a  
163 94 Stockholm  
Tel.: 08 621 34 00  
Fax: 08 621 34 66  
E-mail: info@se.vwr.com

### Switzerland

VWR International AG  
Lerzenstrasse 16/18  
8953 Dietikon  
Tel.: 044 745 13 13  
Fax: 044 745 13 10  
E-mail: info@ch.vwr.com

### UK

VWR International Ltd  
Customer Service Centre  
Hunter Boulevard  
Magna Park  
Lutterworth  
Leicestershire  
LE17 4XN  
Tel.: 0800 22 33 44  
Fax: 01455 55 85 86  
E-mail: uksales@uk.vwr.com